


Applications

Synthesis of carboxylic acids by enzymatic hydrolysis of nitriles.

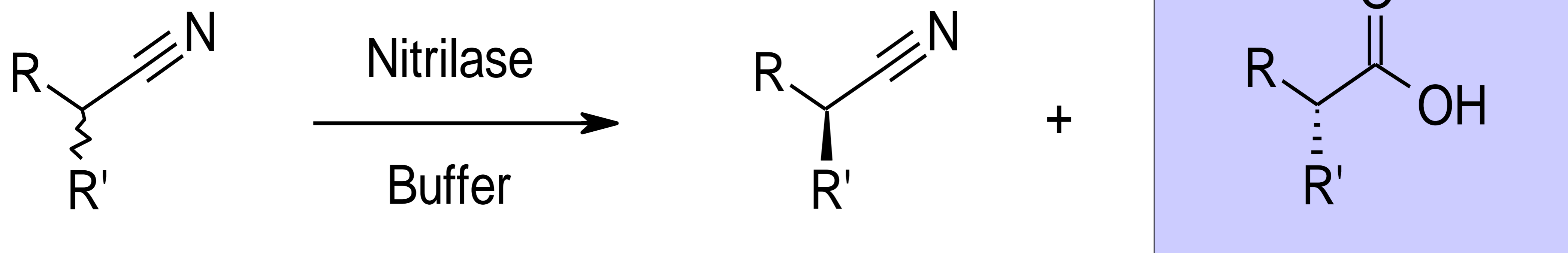

Kit description

The kit contains 96 diverse pre-formulated nitrilase biocatalysts as lyophilised powders in 96-well format, as well as pre-prepared phosphate buffer.

NIT contained in screening kit

	1	2	3	4	5	6	7	8	9	10	11	12
A	101	109	117	125	133	141	149	157	165	173	181	189
B	102	110	118	126	134	142	150	158	166	174	182	190
C	103	111	119	127	135	143	151	159	167	175	183	191
D	104	112	120	128	136	144	152	160	168	176	184	192
E	105	113	121	129	137	145	153	161	169	177	185	193
F	106	114	122	130	138	146	154	162	170	178	186	194
G	107	115	123	131	139	147	155	163	171	179	187	195
H	108	116	124	132	140	148	156	164	172	180	188	196

Contents

Nitrilases	96 enzymes (10 mg each) in 96 well format
DMSO	1 vial (10 mL)
0.1M KH ₂ PO ₄ buffer (pH 7.5)	1 bottle (60 mL)


Screening Procedure

1. Add 500 μ L of the buffer solution to each well*.
2. Add a solution of \sim 1-10 mg substrate in organic solvent (50-150 μ L, depending on solubility) such as DMSO or MTBE to each well.
3. Shake/stir at room temperature (or ideally 30 $^{\circ}$ C). Agitate overnight.
4. Extract product with an organic solvent (MTBE, EtOAc etc.).
5. Analyse sample by chiral GC/HPLC to determine conversion and product ee.

*It is not advisable to keep stock solutions of enzymes, as these will degrade over time. Make each stock solution fresh on the day of use. An adequate supply of nitrilases have been provided for 1 screen.

selectAZyme Offerings

- An ever-expanding biocatalysis team including molecular and microbiologists, enzymologists, bioinformaticians, organic chemists and analysts, all equipped with state-of-the art facilities.
- Expertise in gene identification, expression, fermentation and enzyme production, followed by the efficient use of enzymes to produce complex chiral APIs.
- Enzyme evolution based on computational re-design, semi-rational and random mutagenesis approaches, allowing access to bespoke biocatalysts with enhanced activity, selectivity and process robustness.
- Fully integrated biocatalyst development through screening, (chemo-) enzymatic route definition, process development and scale up (pilot plant facilities available).
- Rapid implementation of enzymatic steps in complex, multi-stage syntheses, leading to significant improvements in production yields and timelines.
- A simple business model that avoids IP issues.

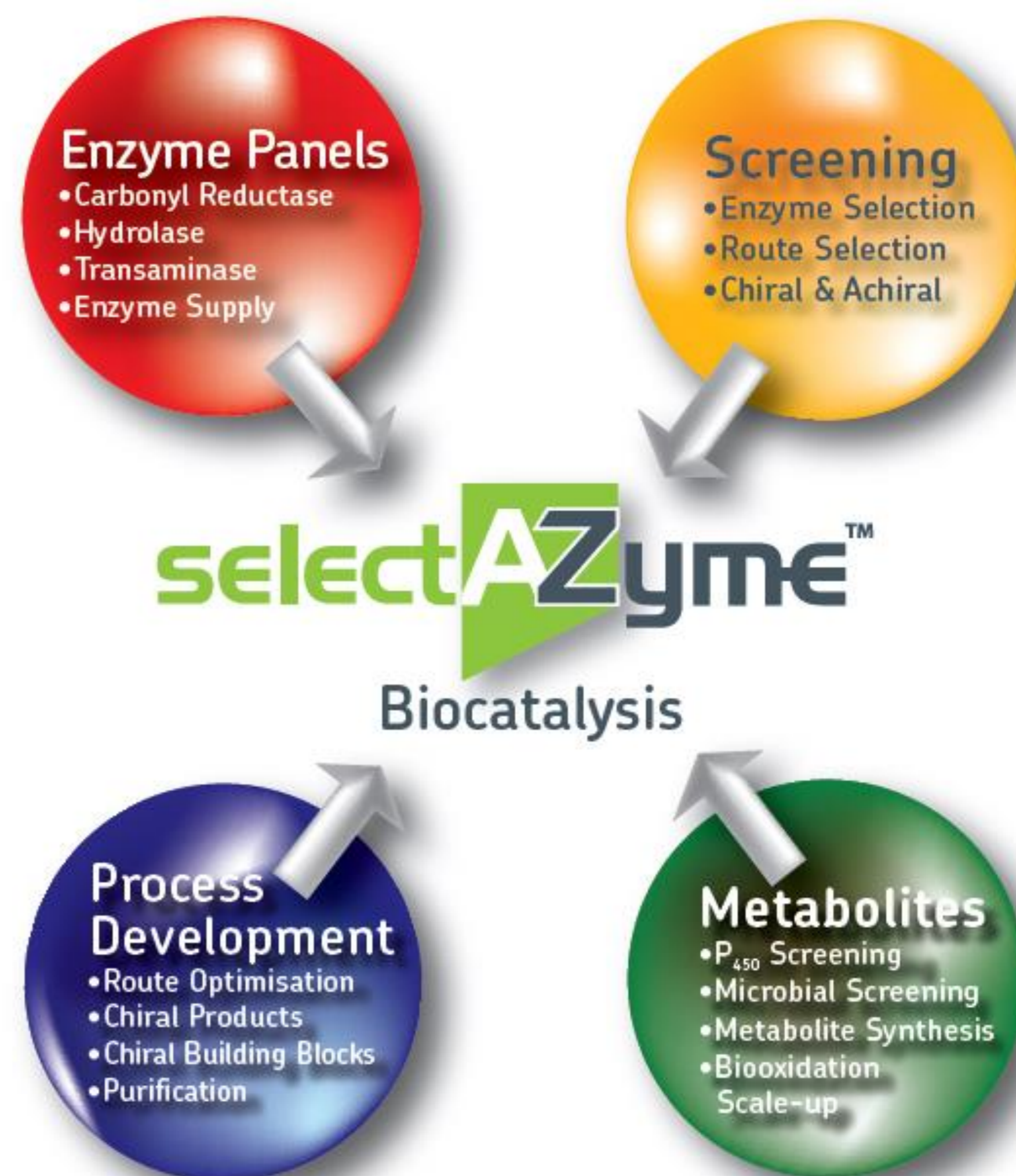

The selectAZyme Range of Enzyme Screening Kits

Our selectAZyme kits include a detailed user guide and come with all buffers, cofactors, recycling systems and reagents necessary to perform screens using standard laboratory equipment.

Carbonyl Reductase (CRED) biocatalysts

96 CRED biocatalysts for the production of chiral alcohols and/or use in cofactor recycling schemes

Aldehyde Reductase (ARED) biocatalysts

16 ARED biocatalysts

Hydrolase biocatalysts

48 commercially available hydrolases for selective acylation of alcohols and amines.

Nitrilase and Nitrile Hydratase (NHase) biocatalysts

9 NHases and 15 nitrilases

Transaminase (TAm) biocatalysts

96 TAm for the production of chiral amines from pro-chiral ketones.

Ene Reductase (ERED) biocatalysts

143 ERED biocatalysts for asymmetric reduction of activated alkenes

P450 Monooxygenase biocatalysts

96 P450 monooxygenase biocatalysts for a huge range of highly selective oxidations

Want Almac to do the screening for you?

- Our experienced biocatalysis team can screen all of our enzymes against your target substrate(s) and simply provide the results.
- Flexible options for subsequent enzyme supply, evolution services, process development and scale up as required.

Technical Contacts:

Prof. Tom Moody, Tel: +44 (0)28 3833 2200 Ext. 5517, E-mail: tom.moody@almacgroup.com.

Dr. Derek Quinn, Tel: +44 (0)28 3833 2200 Ext. 5833, E-mail: derek.quinn@almacgroup.com.

Address: Almac Biocatalysis & Isotope Chemistry Group,

20 Seagoe Industrial Estate, Craigavon BT63 5QD

Web: www.almacgroup.com,

Email: biocatalysis@almacgroup.com